




Calibration Block

The Opentrons Calibration Block is a tool used during Tip Length Calibration and when leveling 8-channel pipettes. It provides a flat surface parallel to the OT-2's deck that all pipettes can reach.

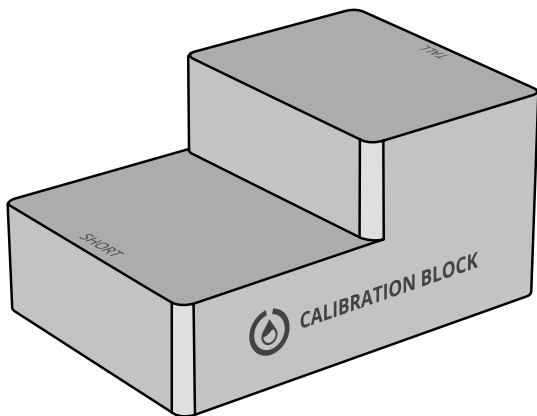

Usage Instructions

To use this Calibration Block with your robot, your Opentrons App and OT-2 must be on software version 4.0 or higher.


If the Opentrons App does not prompt you to use the Calibration Block during Tip Length Calibration, navigate to More > Advanced Settings: Tip Length Calibration Settings. Select “Always use Calibration Block to calibrate.”

For more information about how and when to use the Calibration Block, visit: opentrons.com/support/calibration_block

Support

In case of any issues or questions, contact: support@opentrons.com

Box Contents

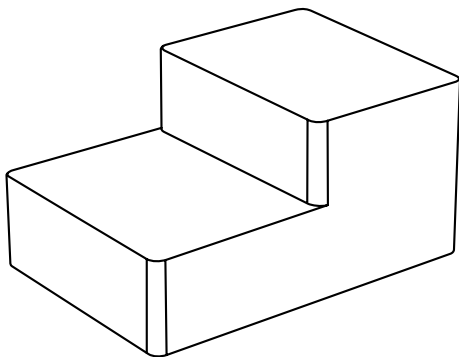

(1) Calibration Block